ITDAF-1

 Tablets must have sufficient mechanical strength to ensure that they will arrive to the end user in the same conditions as they have been produced. One of the most aggressive situations is packing and transport. For this reason, it is convenient to design a test capable to simulate the expected aggressions. For this purpose, two test are developed to measure its parameters **THE FRIABILITY TEST**, and **THE ABRASION TEST**. It consists of a motor that produces a rotatory motion parallel to the working bench; at the end of the motor axle, there is attached a drum, where the samples are introduced, they are weighted before starting the test; when the test is finished, the samples are reweighted again; the weight difference is the weight loss. **FRIABILITY DRUM:** 300 mmØ is made of acrylic plastic, and it is provided with a "Comma" shape internal device, that makes the tablets, do a continuous rolling and falling constant motions, in order to simulate the friability effects, and measure them. **ABRASION DRUM:** 200mm Ø, with 12 blades with 10° angle, make the tablets move up, slide and fall; this motion crumbles them. Metallic painted housing, digital speed control, and timer. Supplied complete with 2 drums (Friability + Abrasion).

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